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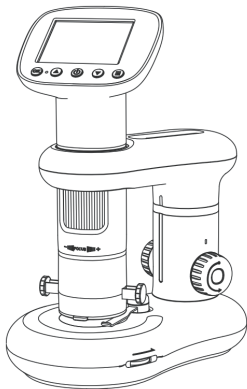
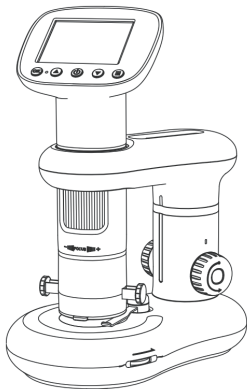
Model: CM3B

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Please read the instructions carefully and keep it for future reference.

# Contents

<b>Read me First</b>	03
<b>1. Installation Instructions</b>	04
Packing list	04
Installation of microscope	04
<b>Parts Description</b>	05
各部件及功能介绍	05
Battery installation	06
<b>3. Detailed Operation</b>	07
How to focus?	07
How to work with PC?	08
App download	08
Operating steps	09
USB function description	10
<b>4. Interface Introduction</b>	11
Photo mode	11
Video mode	11
Playback mode	12
Menu interface	13
Function introduction	14
Usage Environment Description	
<b>5. Product Parameters</b>	16
<b>6. 注意事项</b>	17

# Read me First

- Please read this user manual carefully and keep it for future reference.
- Pay attention to protect the lens part. Close the lens lid after using to avoid entering dust or strange objects.
- The microscope can be used in hand or inserted as a desktop, with attention to avoid damage to the components.
- When placing items, place them directly below the center of the lens as much as possible. Avoid not seeing objects clearly and bringing about a bad experience.
- Please store the microscope in a cool and dry place and avoid direct sunlight.
- Please avoid storing the microscope in places where there is risk of falling, which may cause the internal lenses to deviate or damage other components, as well as other irreparable damage.

---

---

① **Note: The plan used in this manual is for reference only, please refer to the actual plan.**

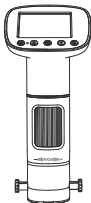
---

---


# 1 .Installation Instructions

- Packing list

Microscope

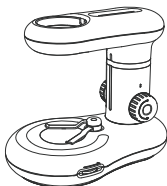Base

- Installation of microscope

1. Take off the microscope lens cover and transparent compartment .


2. Hold the base with one hand and insert the microscope vertically into the installation hole.

3. Install the transparent compartment, Complete

## 2 .Parts Description
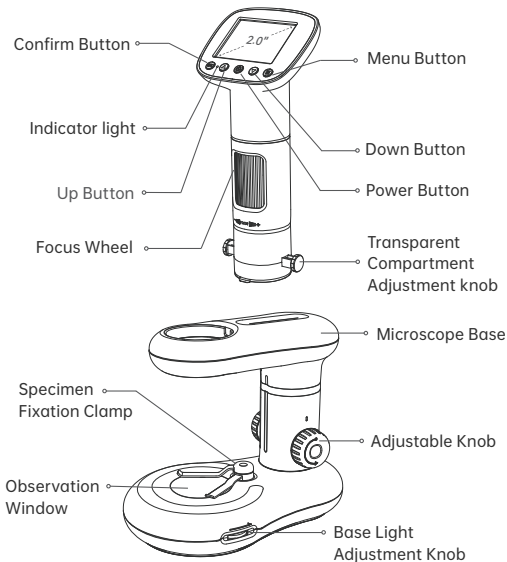
### • Parts and functions

### **OK/Confirm Button**

In the Mode interface, it is the Photo/Video/Play button. On the Menu interface, it is the Confirm button

### **Menu Button**

Used to indicate charging, computer connection and working status

### **Indicator light**

Used to indicate charging, computer connection

### **Power Button**

Long press 3 seconds to turn on/off; Press to adjust the brightness

### **Up Button**

Switch up

### **Down Button**

Switch down

### **Specimen Fixation Clamp**

A clamp for fixing specimens or observing targets

### **Observation Window**

Render the bottom of the object for observation

### **Adjustable Knob**

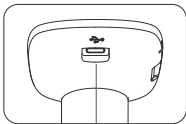
Adjustable Movement of Microscope for Up and Down

### **Base Light Adjustment Knob**

Rotate the knob to adjust the brightness of the base light.

## • Parts and functions

### Type-C Interface

To charge or connect with computer.

### Reset Button

When the microscope is abnormal, poke this hole with a fine pointed object to force the power off.

### TF card slot

Support maximum 32G memory card (Note: please use C10 high-speed card)

You can charge the microscope with your daily mobile charger, computer, or power bank.

Please do not charge it for more than 12 hours to avoid affecting the battery life.

---

**① Note : During the charging process, if the outer shell of the microscope has slight heating, it is a normal phenomenon, Please use it with confidence.**

---

## • Battery installation

**Step 1.** Pry open the battery cover of the base by buckling the battery cover recess

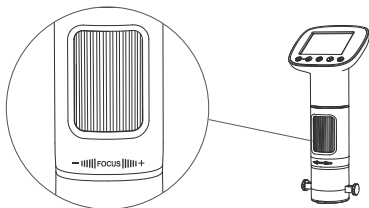
**Step 2.** Install the 7# 1.5V battery according to the direction indicated inside (please pay attention to distinguish the positive and negative poles when installing the battery, the non-rechargeable battery can not be charged )

**Step 3.** Close the battery cover.

### 3 .Detailed Operation

- ① Before using the microscope, please make sure to hold it steady with your hands or place it on the desk for use.

#### • How to focus?

-  Counter-Clockwise Rotation: multiple magnification
-  Clockwise Rotation: multiple reduction
- ( **Note:** In case of any other problems, the reset hole will restore the factory setting. At the same distance, the rotating zoom cylinder has two magnification. To adjust more multiples, it can be adjusted up and down through the microscope base bracket. )

## • How to work with PC?

### 1.App download

◉ **Note: The following 3 methods can be downloaded, please choose the appropriate version according to your computer system.**

① For Windows 7, Windows 8.1, Windows 10, Windows 11 and Mac , download the software from the website:  
<http://soft.hvscam.com>

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
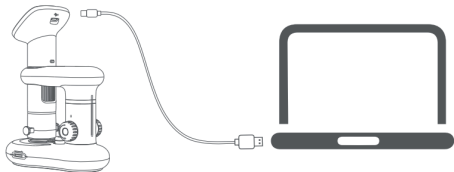
③ For Mac OS X 12 or higher system, search "hiView" in the APP Store to download it and you can also be viewed in Photo Booth(“”).

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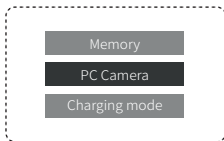
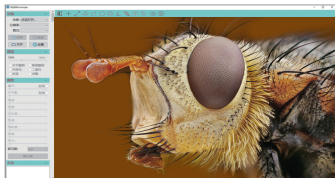
### 3.USB function description

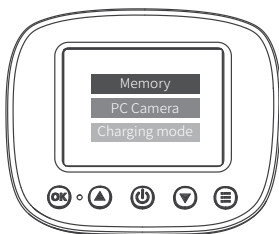When connecting the device to a computer through a USB cable, there are 3 modes of Memory/PC Camera/Charging mode available to the user.

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③ **Charging mode:** When the power of the microscope is insufficient, you can use the microscope while charging, connect the microscope and the computer with the Type-c data cable, start the microscope, select the charging mode, you can enter the charging state.

---

④ **Note:** You can also charge the microscope with your daily phone charger, computer or power bank.

---


## 4.Interface Introduction

- ⓘ (Note:Before using the microscope ,please insert the TF card into the microscope to ensure better storage of Photo/Video file)


### • Photo mode

- If the icon in the upper left corner of the microscope is "📷", the current mode is Photo mode. Long press the Menu button on the microscope under the startup state to switch to the Photo mode.

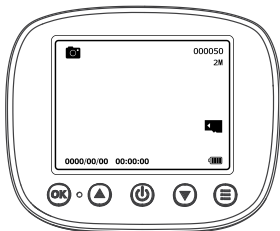
### • Video mode

- If the icon in the upper left corner of the microscope is "👤", the current mode is Video mode. Long press the Menu button on the microscope to switch to the Video mode under the startup state.
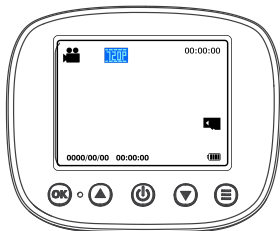
### • Playback mode

- If the icon in the upper left corner of the microscope is "▶", the current mode is Playback mode. Long press the Menu button on the microscope to switch to Playback mode under the startup state.

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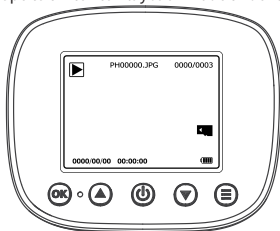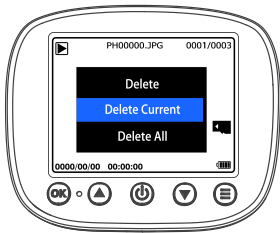
## ● Function introduction

- ◎ **Photo resolution:**有2M,3M,5M三种格式可选择。
- ◎ **Video resolution:** there are 1080P, 720P, 2 kinds of formats can be selected.
- ◎ **时间标志:** select “On”,the lower right corner of the photo/video file will display the date and time logo; select “off”,the photo/video file does not display the date and time logo.
- ◎ **Language settings:** 英文, 简体中文, 繁体中文, 日语, 韩语, 俄语, 法语, 德语, 葡萄牙语, 波兰语, 荷兰语, 希伯来语, 意大利语, 西班牙语)。
- ◎ **屏幕亮度:**有1-8共8个亮度挡位可供选择。
- ◎ **Automatic shut-down:**select the auto shutdown time, the microscope will be automatically shut down when no one is operating; no auto shutdown, please select "Off".
- ◎ **Screen Protector:** select the appropriate time to make the device automatically turn off the screen when no one is operating.
- ◎ **Format:**This operation will delete all data in TF card. (注:删除后将无法恢复, 请谨慎操作)
- ◎ **Date/Time:**You can set the date / time parameter of microscope display.
- ◎ **Restore the factory settings:** 所有设置都会恢复到出厂默认参数。
- ◎ **Version:** You can view the current firmware version.
- ◎ **Delete:** 在回放模式查看单个图片或视频时,短按菜单键可将相册内文件进行删除,可选择删除当前或全部删除。

## • Usage Environment Description

**ⓘ (Note: The capacity of the battery may be reduced to some extent in low temperature conditions.)**

Please use in an environment of -15°C~55°C.

Avoid wetting the microscope with water, drinks, and corrosive liquid.

Avoid keeping the microscope close to heat sources, open flames, and flammable or explosive gases (or liquids).

Avoid dust entering the lens and interior of the microscope.

Avoid impacts and strong vibration on the microscope.

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